



GlobusWorld 2024

Automating Instrument Data at Scale

Rachana Ananthakrishnan
Vas Vasiliadis






Data processing pipeline pattern



Data source

At the minimum, make data available to collaborators

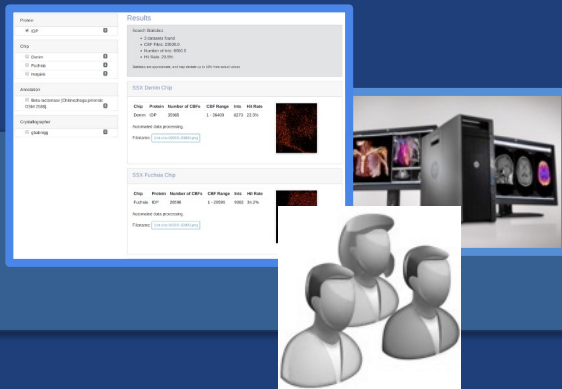


Processing and storage



Set permissions for access

Access by collaborators

g Data processing pipeline pattern

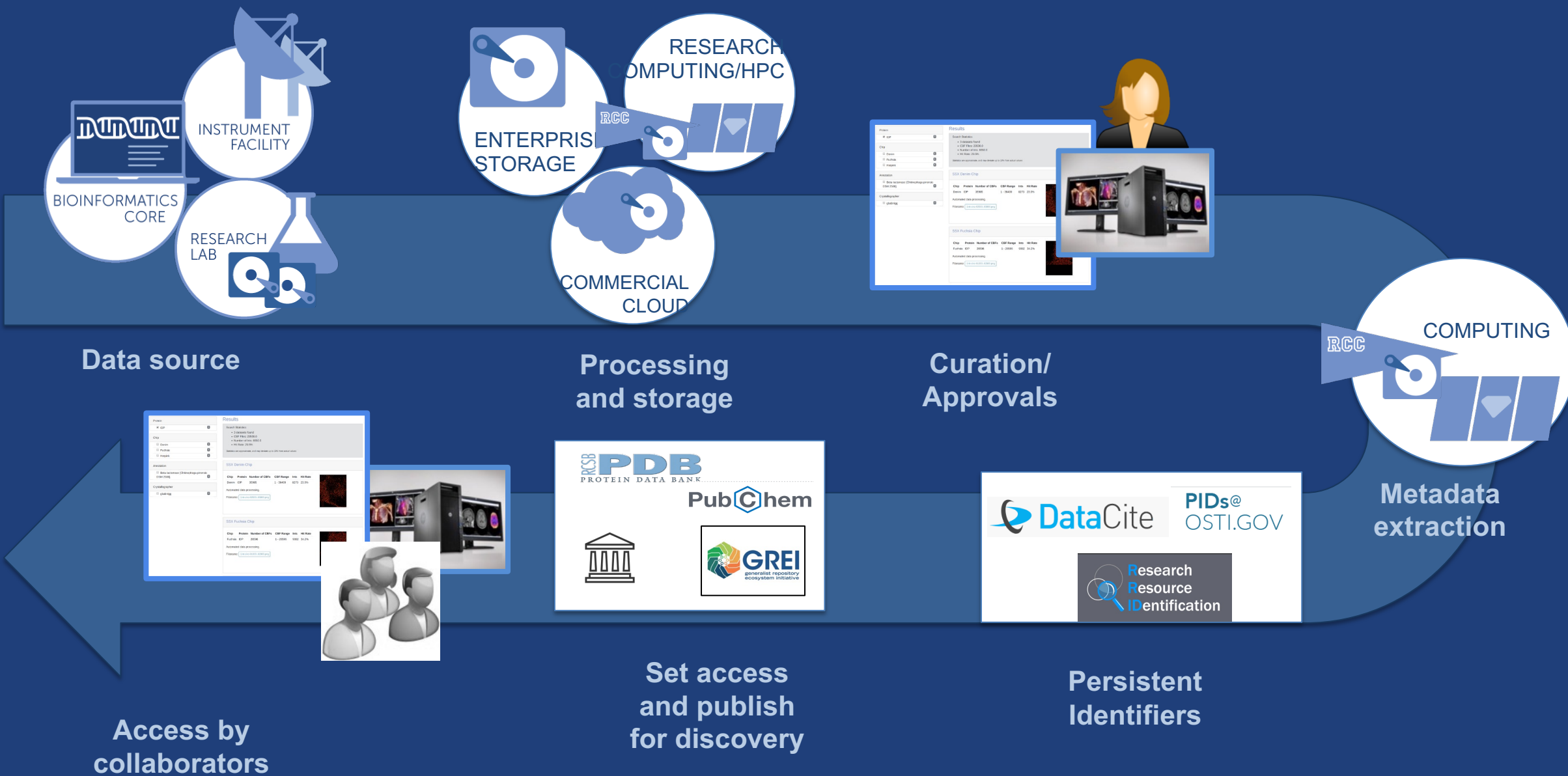

Requirements for such pipeline

- **Reliable, near-real time data access**
- **Uniformed policy for data access, based on local policy**
- **Delegation of data access management to PI**
- **Ability to compute on data across storage classes**
- **Apply best practices with data processing pipeline**
- **Support for data organization to facilitate FAIR data**
- **...**


Automation to scale



How do we use the Globus platform for full automation?



Data processing pipeline pattern

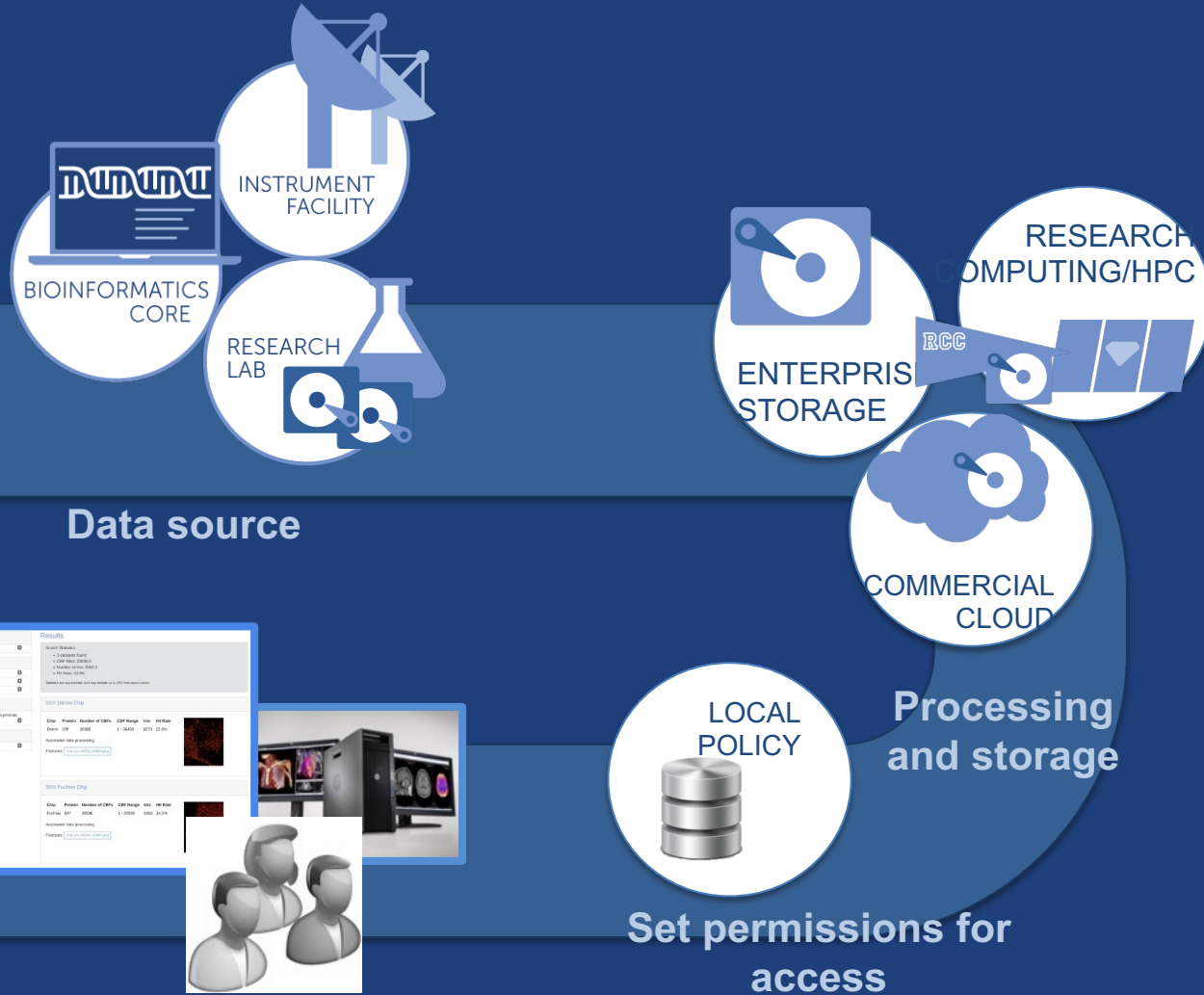

1. Preparing acquisition machine for data access
2. Configuring data collections for automated data transfer
3. Credentials for automation
4. Create and deploy flow for automation
5. Create trigger scripts



1. Preparing acquisition machine for data access



Install Globus Connect

- **For Linux machines, install and use Globus Connect Server**
- **Acquisition machines often are Windows machine**
 - **Globus Connect Personal deployment**
 - Installed as a local user account
 - PI login vs GCP always running on a shared account
 - Outbound connections only



2. Configuring data collections for automated data transfer

Using guest collections at instrument facilities

- **Create guest collections on all storage systems for automation**
 - Acquisition machine, storage mounted on compute etc.
- **Creation of guest collection cannot generally be automated ***
- **Permission and role management on guest collections can be fully automated**



Pattern 1: Guest collection use

- **Create a guest collection at top level directory**
 - This is done by a user who has local account
- **For each experiment/project/modality**
 - Create a folder
 - Set permissions PI/collaborators to read data from the folder
- **Can automate permission management by using local policy store**

Pattern 2: Guest collection use

- **Create a guest collection for each experiment/project/modality**
 - Grant PI role to manage access to the guest collection (Access Manager role)
 - Set permissions collaborators to read data from the folder
- **Can automate role and permission management by using local policy store**



3. Credentials for automation


Managing service accounts/app credentials



- **Application Identity:**
appclientid@clients.auth.globus.org
- **These are confidential apps with client id and secret**
- **Ensure application is on a secure device**
- **Set up policy for rotation of secret**
- **Assign project admins to manage the registration**

Registering a service account

- **Webapp - Settings**
 - app.globus.org/settings/developers

Register an...

App 

 Register a service account or application credential for automation 

Applications that authenticate and act as the application itself. These applications are used for automation and as service or community accounts, and do NOT act on behalf of other users.



Get app credentials at app.globus.org/settings/developers

Application Registration

App Registration

Project Name: Client Type Tests

App Name*
A human-friendly name to help you keep track of this app.

Native App Will be used by a native application

Redirects: `https://example.com/authcallback`
One URL per line – must be HTTPS

Required Identity Require that the user has linked an identity from a specific Identity Provider, but does not require that the user authenticate with that identity each time.

Pre-select Identity Provider Pre-select a specific Identity Provider on login page

Use effective identity (ID token + userinfo) Identity returned in userinfo is the primary identity unless a required identity is selected above. If unchecked, the identity returned in userinfo is the most recently authenticated identity. (This should usually be checked for most applications.)

Privacy Policy URL: `https://example.com/privacy`

Terms & Conditions URL: `https://example.com/terms-and-conditions`

[Register App](#) [Cancel](#)

[Edit App Details](#)

[+ Add Client Secret](#)

[Delete App Registration](#)

This secret will only be shown once, please copy and store it securely. If lost, Globus will not be able to recover it, and you will need to register a new secret.

Secret

`z8pQAik84azJXLWys/vZWYX0jqXGOxY3+vRsLuM7lbg=`

[Done](#)



Get app credentials at

app.globus.org/settings/developers

The screenshot shows the 'Settings' page in the Globus application. The 'Developers' tab is selected in the top navigation bar. The main content area is titled 'Register an...' and lists several options for registering different types of applications, each with a brief description and an icon.

- App**: Register a service account or application credential for automation. Applications that authenticate and act as the application itself. These applications are used for automation and as service or community accounts, and do NOT act as users for other users.
- Portal**: Register a portal, science gateway, or other application you have. Applications that are OAuth confidential clients, and can manage a client secret on behalf of users.
- Thick client**: Register a thick client or script that will be installed and run by users on their devices. Applications that are OAuth public clients or native applications that are distributed to users, and thus cannot manage a client secret.
- Service API**: Register a service API that will use Globus for authentication. A RESTful API (OAuth resource server) that uses Globus Auth tokens for authentication. Users accessing the service login via Globus and consent for the client to use the service.
- Advanced registration**: Register any client with Globus Auth - an application (confidential or public), a service account, or service API.

The screenshot shows the details for a registered application named 'Django ESGF2 Portal'. The page includes a sidebar with navigation options (Apps, Policies, Roles) and a main content area with the following details:


- Client UUID**: b19fbd77-a940-47bb-98d3-0fdec16c0bc6
- Identity**: b19fbd77-a940-47bb-98d3-0fdec16c0bc6@clients.auth.globus.org
- App Type**: Globus Connect Server
- Client Secrets**: A table listing secrets for the application.

Client Secret	Created	UUID
Search and auth clients for Globus SDK	Created 234 days ago	UUID: bc7ad45d-4bdd-4854-8b5d-52b4b38583b0

On the right side of the page, there are three action buttons: 'Edit App Details', 'Add Client Secret', and 'Delete App Registration'.



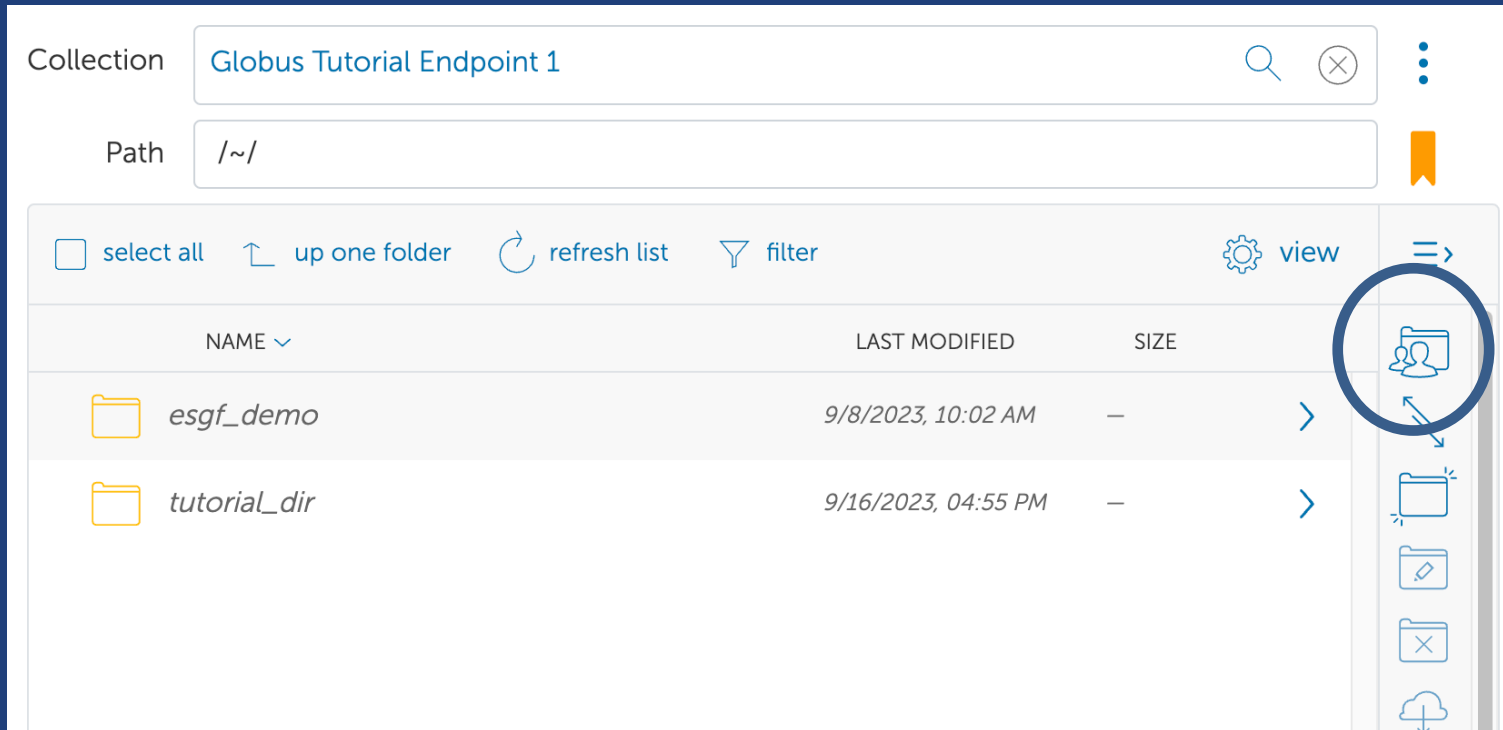


Set permission for the service account






- Use the web app to create guest collection on guest collection you created









Collection: Globus Tutorial Endpoint 1

Path: /~/

select all  up one folder  refresh list  filter  view 

NAME 	LAST MODIFIED	SIZE	
 <i>esgf_demo</i>	9/8/2023, 10:02 AM	–	
 <i>tutorial_dir</i>	9/16/2023, 04:55 PM	–	

Vertical toolbar icons (from top to bottom):
 (circled in blue)








Set permission for the service account



Overview Permissions Roles

Assigned Roles

[Assign New Role](#)

	Rachana Ananthakrishnan	ranantha@uchicago.edu	Administrator
--	-------------------------	-----------------------	---------------

Assign To:

Role

- Administrator
modify endpoint definition, delete the endpoint, manage roles, perform file system operations and transfers, and all capabilities of the Access Manager and Activity Manager roles
- Access Manager
view, add, and delete all access rules on the endpoint; implicitly gives read/write access to the root of the endpoint
- Activity Manager
view and control tasks and other endpoint activity
- Activity Monitor
view tasks and other activity to or from the endpoint

[Add Role](#) [Cancel](#)



4. Create and deploy flow for automation



Create a flow for your use case

- **Start with flow definitions that are published:**
 - github.com/globus/globus-flows-trigger-examples
 - docs.globus.org/api/flows/authoring-flows/examples/
- **Manage flow definitions in a version controlled system**
- **Validation tools**
 - Flows IDE: <https://globus.github.io/flows-ide/>
 - Globus CLI: *globus flows validate*



Flow definition

```
"StartAt": "TransferFiles",
"States": {
  "TransferFiles": {
    "Comment": "Transfer to a guest collection",
    "Type": "Action",
    "ActionUrl": "https://actions.automate.globus.org/transfer/transfer"
    "Parameters": {
      "source_endpoint_id.$": "$.input.source.id",
      "destination_endpoint_id.$": "$.input.destination.id",
      "transfer_items": [
        {
          "source_path.$": "$.input.source.path",
          "destination_path.$": "$.input.destination.path",
          "recursive.$": "$.input.recursive_tx"
        }
      ]
    },
    "ResultPath": "$.TransferFiles",
    "WaitTime": 60,
    "Next": "SetPermission"
  },
  "SetPermission": {
    "End": True
  }
}
```

Action

Action Provider URL

Action inputs

Timeout (seconds)

Next state



Choice in flows

```
"DetermineOutcome": {  
  "Type": "Choice",  
  "Choices": [  
    {  
      'Variable': "$.curator_decision",  
      'StringEquals': "approve",  
      'Next': "SetPermissionForAccess"  
    },  
    {  
      'Variable': "$.curator_decision",  
      'StringEquals': "approveEmbargo",  
      'Next': "SetPermissionForEmbargo"  
    }  
  ],  
  "Default": "SubmissionRejected"  
},
```

Action Type

Evaluate and
next state

Evaluate and
next state



Run context

- **Flow and user properties as read-only values**
- **Available in `$_context`**
 - Flow id
 - Run id
 - All identities of the user invoking the flow
 - Email id of the user invoking the flow
 - Token information (issued time, and expiration information)






Permissions to run the flow

Assign To

- User
- Group
- All Logged In Users
- Public (anonymous users)

Role

-  Administered By
can start this flow, view this flow and associated activity, and modify this flow
-  Runnable By
can start this flow and view associated activity
-  Visible To
can view this flow and associated activity


Add Role

Cancel


**Set
permission
for the service
account to
run the flow**



Triggering flows on instruments ...and other resources

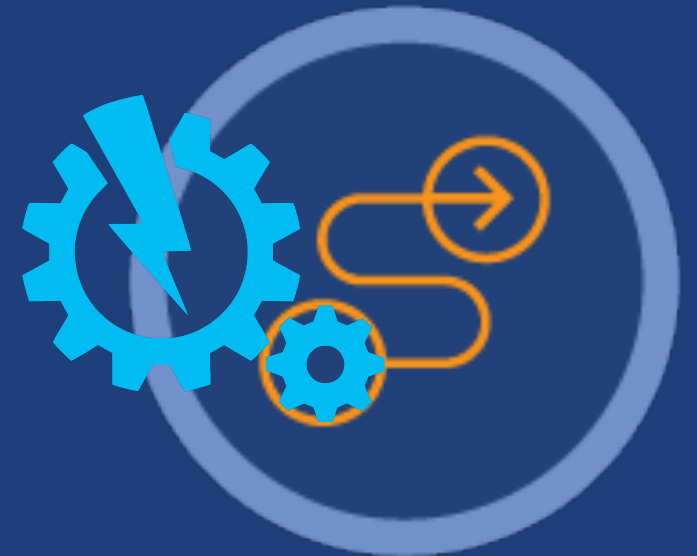

 A simple, and very common, use case

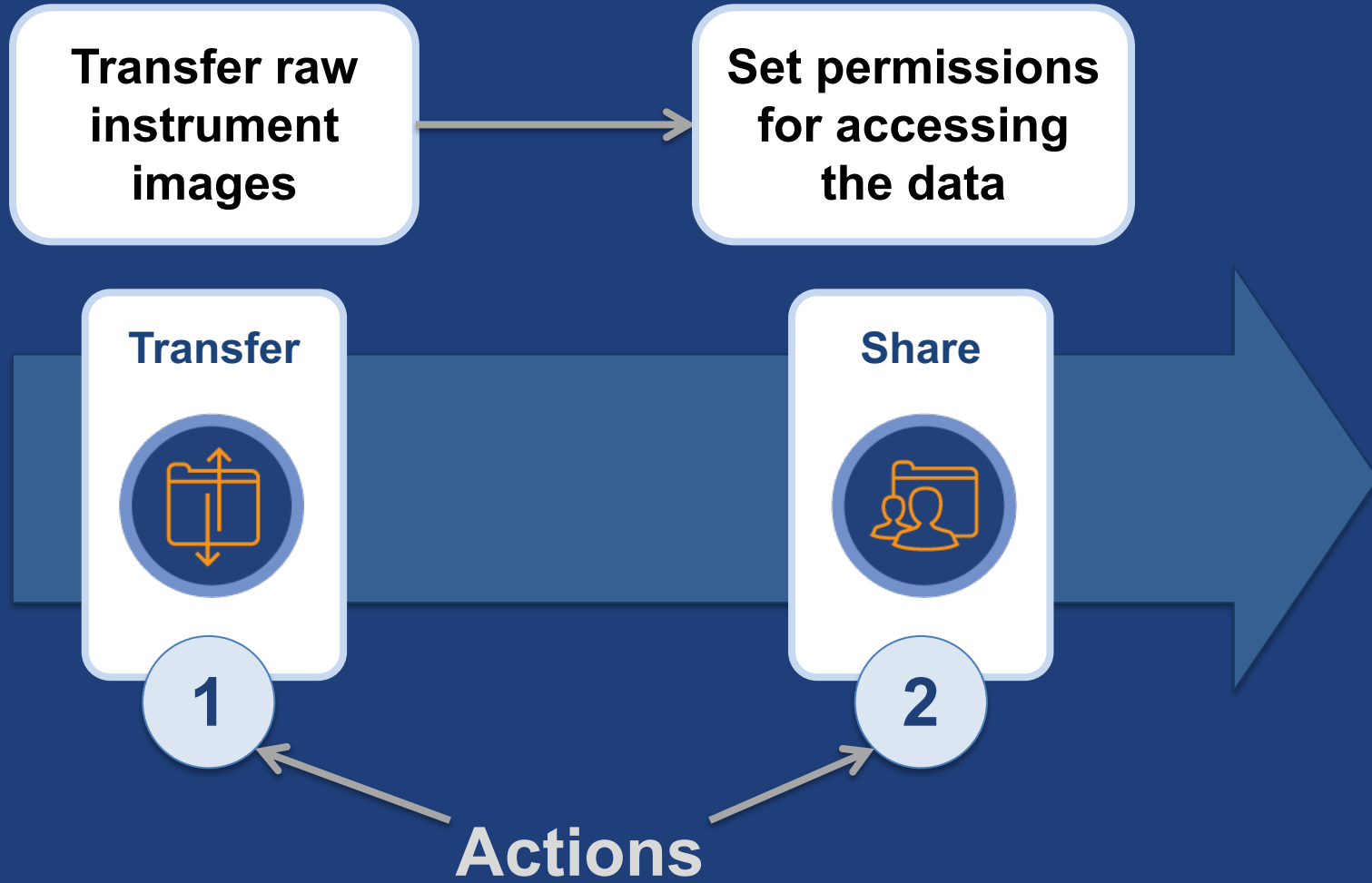

Creating and triggering runs of this flow



1. **Log into the Globus CLI**
2. **Create (deploy) the transfer-and-share flow**
3. **Edit the monitor (trigger) script**
4. **Ensure GCP is running on the instrument**
5. **Run the monitor script**
6. **Trigger the flow**

github.com/globus/globus-flows-trigger-examples



Create the flow



```
$ globus login
$ cd ~/globus-flows-trigger-examples/transfer_share
$ globus flows create FLOW_NAME \
> definition.json --input-schema schema.json
$ ~/globusconnectpersonal-3.2.5/globusconnectpersonal -start &
```

- Success returns the flow ID
- Inspect the flow using the web app

Run the monitor script and trigger the flow



```
$ source ~/.trigger/bin/activate
$ cd ~/globus-flows-trigger-examples
$ ./trigger_transfer_share_flow.py \
> --watchdir /home/dev3/images \
> --patterns .done
$ cp ~/test-data ~/images
$ touch ~/images/i.am.done
```


Directory to monitor for file creation

Flow is triggered when a new filename matches this expression

Simulate instrument data creation

Trigger the flow

Adding computation to our instrument flow

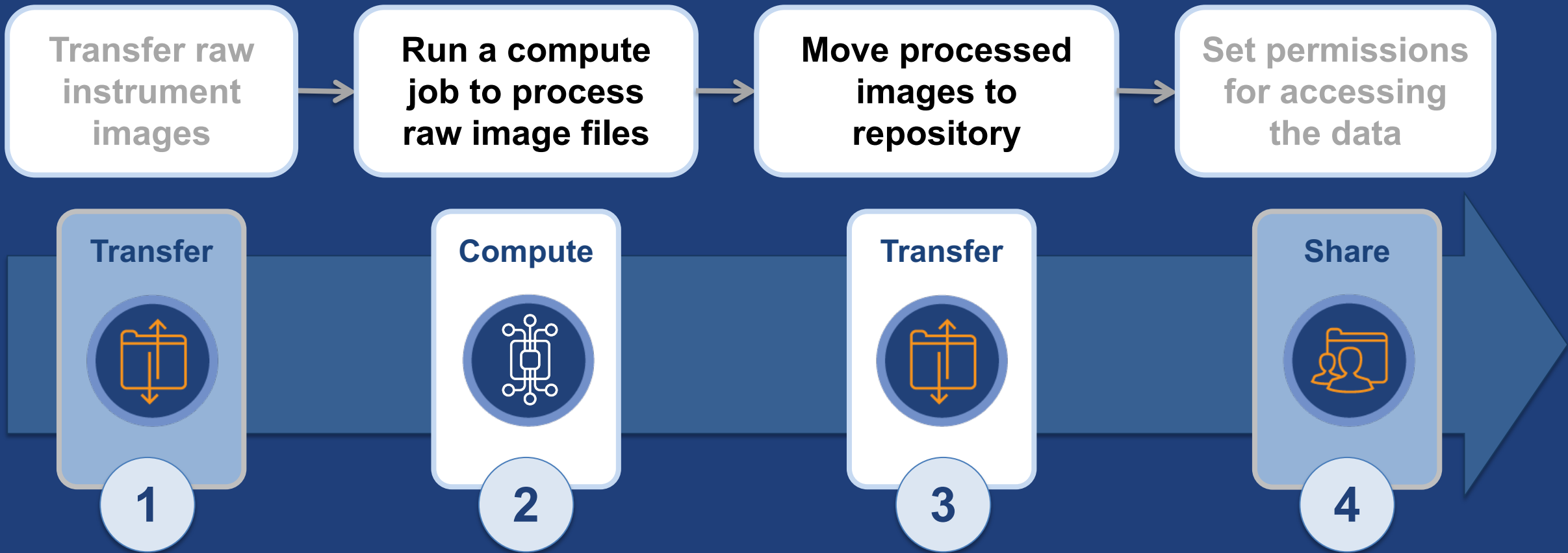

Our instrument research environment

Registered
Compute
Function

Sharing Repository
(ALCF Eagle)



access
result files

Storage
Endpoint

HPC System
(UChicago RCC Midway3)



Compute
Endpoint


```
def process_images(input_path=None, result_path=None):  
    import os  
    import glob  
    from PIL import Image  
  
    files = (file for file in glob.glob('*.*')  
            if os.path.isfile(os.path.  
                if not os.path.exists(result_path):  
                    os.makedirs(result_path)  
  
    for file in files:  
        image = Image.open(file)  
  
        # Generate thumbnail  
        image.thumbnail((200, 200))  
  
        # Save thumbnail image  
        image.save(f'{result_path}/thumb_{file}')  
    
```



invoke image
processing function



"The Instrument"
(just a VM in the cloud :-)

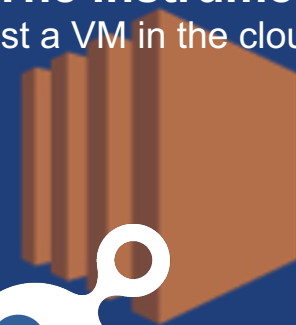

Monitor
script



trigger
flow run

Storage
Endpoint
(using GCP)

transfer
raw files



transfer
result files




set
permissions



Transfer
Service



transfer control

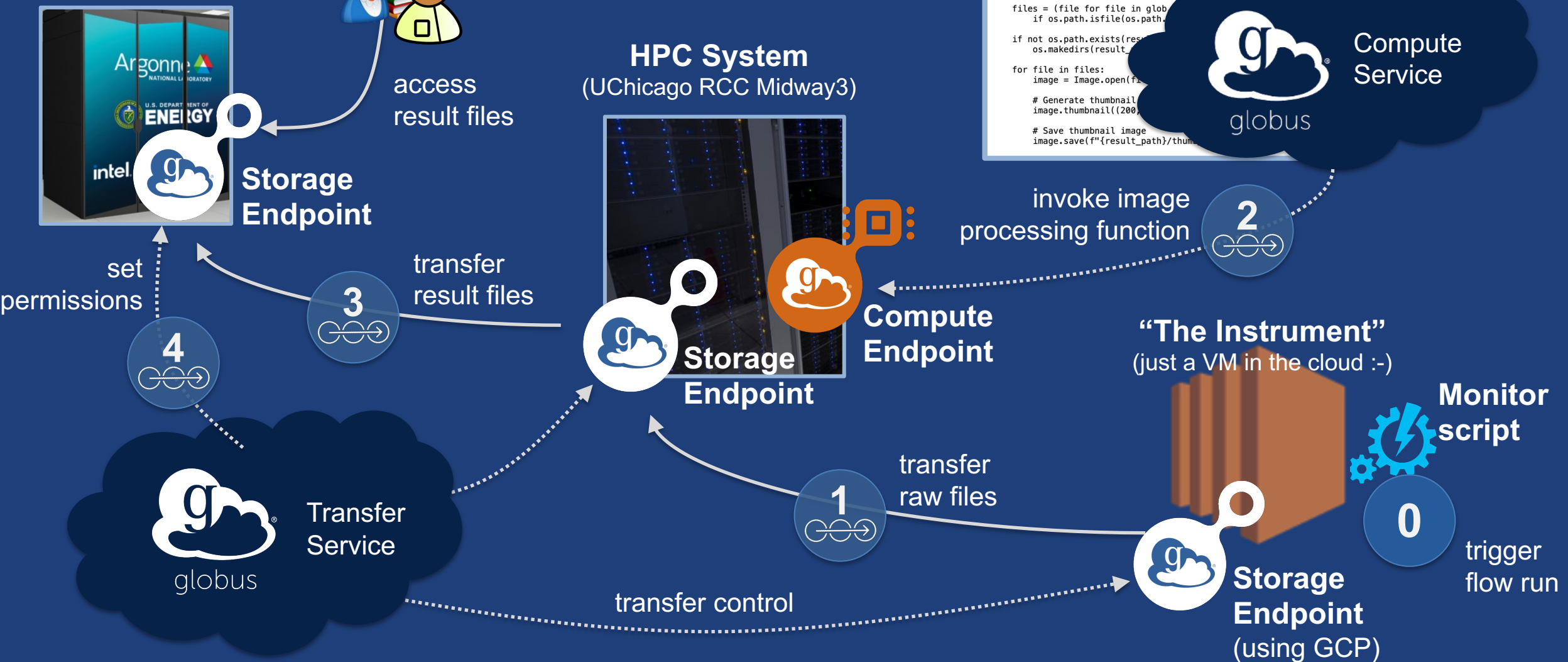



Key consideration: Be identity aware



- **What identity is the flow running as?**
- **Does identity have access to target resources?**
 - Collections (ideally, guest collections)
 - Compute endpoint
 - Compute function
- **Does identity have the required role?**
 - Access Manager, if granting/revoking permissions



Making Data Findable with Globus Search



Data description and discovery

- **Metadata store with fine-grained visibility controls**
- **Schema agnostic dynamic schemas**
- **Simple search using URL query parameters**
- **Complex search using search request document**



docs.globus.org/api/search



Data ingest with Globus Search

POST /index/{index_id}/ingest'

```
{
  "ingest_type": "GMetaList",
  "ingest_data": {
    "gmeta": [
      {
        "id": "filetype",
        "subject": "https://search.api.globus.org/abc.txt",
        "visible_to": ["public"],
        "content": {
          "metadata-schema/file#type": "file"
        }
      },
      ...
    ]
  }
}
```



- Bulk create and update
- Task model for ingest at scale



Data ingest with Globus Search

POST /index/{index_id}/ingest'

```
{
  "ingest_type": "GMetaList",
  "ingest_data": {
    "gmeta": [
      {
        "id": "weight",
        "subject": "https://search.api.globus.org/abc.txt",
        "visible_to": ["urn:globus:auth:identity:46bd0f56-
          e24f-11e5-a510-131bef46955c"],
        "content": {
          "metadata-schema/file#size": "37.6",
          "metadata-schema/file#size_human": "<501b"
        }
      },
      ...
    ]
  }
}
```



Visibility limited to Globus Auth identity

- Single user
- Globus Group
- Registered client application

Data discovery with Globus Search

GET /index/{index_id}/search?q=type%3Ahdf5

Simple query

```
{
  "@datatype": "GSearchResult",
  "@version": "2017-09-01",
  "count": 1,
  "gmeta": [
    {
      "@datatype": "GMetaResult",
      "@version": "2019-08-27",
      "entries": [
        { ... }
      ],
      "subject": "https://..."
    }
  ],
  "offset": 0,
  "total": 1
}
```




Data discovery with Globus Search

POST `/index/{index_id}/search`

```
{
  "filters": [
    {
      "type": "range",
      "field_name": "pubdate",
      "values": [
        {
          "from": "*",
          "to": "2020-12-31"
        }
      ]
    }
  ],
  "facets": [
    {
      "name": "Publication Date",
      "field_name": "pubdate",
      ...
    }
  ]
}
```

Complex query

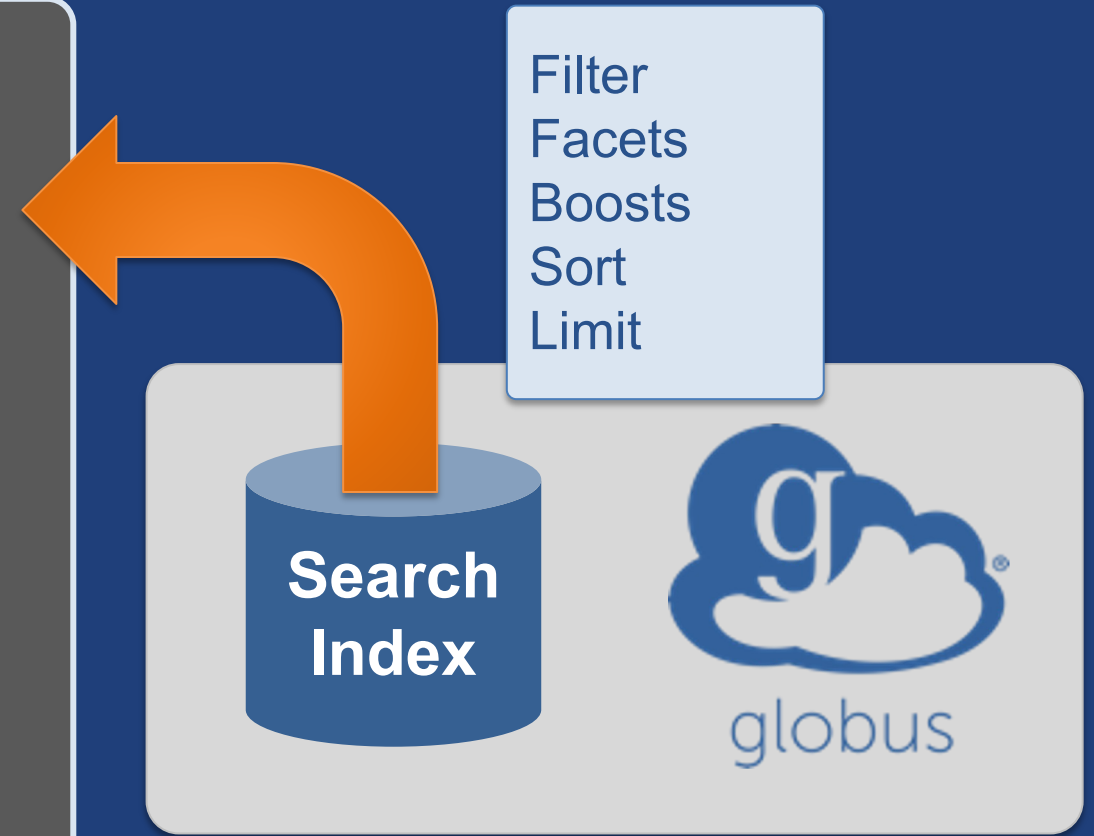



Making our instrument data FAIR

Transfer raw instrument images

Run a compute job to process raw image files

Move processed images to repository

Transfer



Compute



Transfer



Share



Search



Search



Set permissions for accessing the data

Ingest protected metadata, searchable by one group

Ingest open metadata, searchable by all

End-to-end automation in practice: XPCS

Data capture



Globus
Flows



Transfer



Transfer
HDF5 files

Transfer



Transfer
IMM

Compute



Run Corr



FAIR data, ready for discovery!

Protein

IDP

Chip

Denim

Fuchsia

Hotpink

Annotation

Beta-lactamase [Chitinophaga pirinis DSM 2588]

Crystallographer

gbabnigg

Results

Search Statistics

- 3 datasets found
- CBF Files: 23500.0
- Number of Ints: 6650.0
- Hit Rate: 29.5%

Statistics are approximate, and may deviate up to 10% from actual values

SSX Denim Chip

Chip	Protein	Number of CBFs	CBF Range	Ints	Hit Rate
Denim	IDP	35985	1 - 36400	8273	23.0%

Automated data processing.

Filename: `3ire-sinc-02000_02000.png`

SSX Fuchsia Chip

Chip	Protein	Number of CBFs	CBF Range	Ints	Hit Rate
Fuchsia	IDP	26596	1 - 26596	9082	34.2%

Automated data processing.

Filename: `3ire-sinc-01000_02000.png`

Search



Ingest to
index

Share



Set access
controls

Transfer




Move results
to repo

Compute




Gather
metadata

Compute

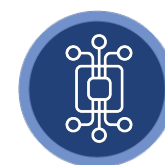

Plot results

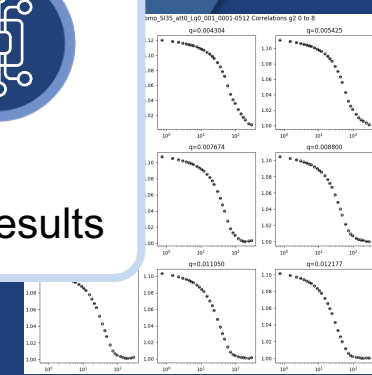



Take a look...

acdc.alcf.anl.gov

Extending the ecosystem: Action Providers

- **Action Provider is a service endpoint**

- Run
- Status
- Cancel
- Release
- Resume



docs.globus.org/api/flows/hosted-action-providers

- **Action Provider Toolkit**

action-provider-tools.readthedocs.io



Custom developed